

AFTER REACTION IS FINISHED

- TIP UPPER LAYER FROM SIDE-ARM BACK INTO RB FLASK CAREFULLY, LEAVING HOH IN SIDE-ARM.
- PIPET ESTER INTO CENTRIFUGE TUBE CONTAINING HOH. ADD TO THIS CONTENTS OF SIDE-ARM.
- MIX CONTENTS WELL, REMOVE HOH LAYER, EXTRACT ORGANIC LAYER WITH AQUEOUS BICARBONATE, THEN SAT'D NaCl.
- DRY ORG LAYER WITH CaCl₂.
- DISTILL TO PURIFY ESTER.
- WEIGH, CALC YIELD, TAKE IR SPECTRUM.